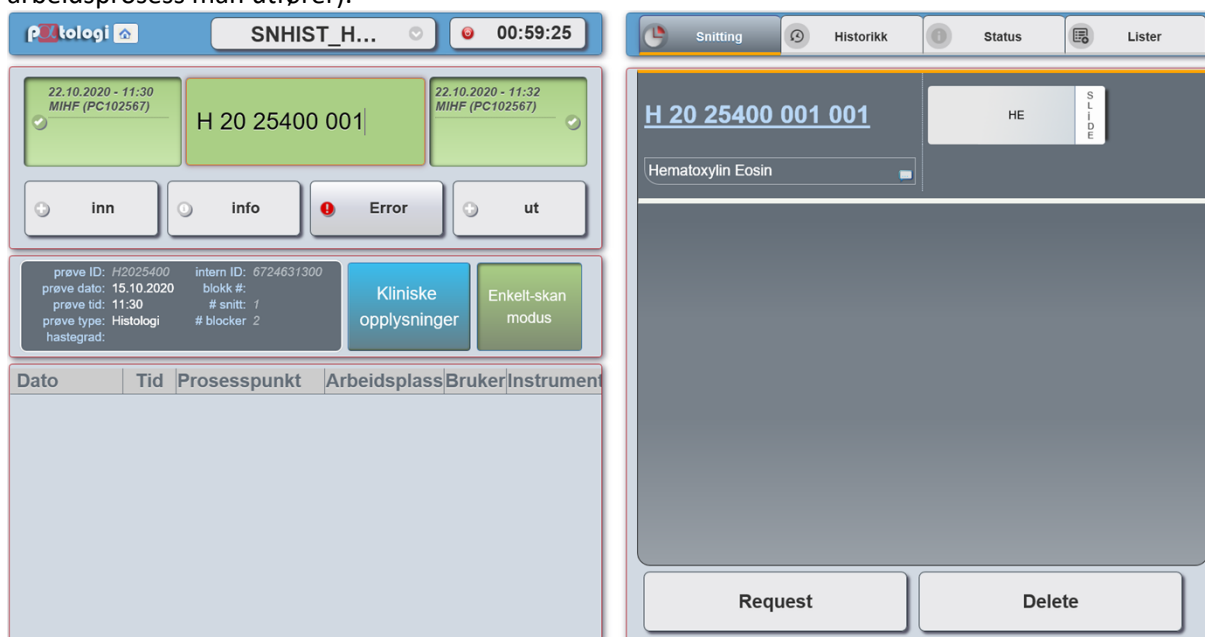
 Helse Bergen HF/Laboratorieklinikken Kategori: []	<h2>Unilab - UniApp / prosesser (PAT)</h2>
	Brukerveiledning
Dok. eier: <a href="#">Johanne Lind Aasen</a>	Dok. ansvarlig: <a href="#">Cecilie Norheim</a>
Organisasjon: Helse Bergen HF/Laboratorieklinikken	Dok.ID: D65096
Kategori: []	Versjon: 2.01/06.02.2024 Gyldig til: 06.02.2025

## 1 UniApp / prosesser

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The image shows two screenshots of the UniApp interface. The left screenshot displays a dialog for a process point with a timer at 00:59:25. It includes fields for 'inn', 'info', 'Error', and 'ut', and a section for 'Prøve informasjon' with fields for preve ID, preve dato, preve tid, preve type, hastegrad, intern ID, blokk #, # snitt, and # blocker. The right screenshot shows a detailed view of a process point with ID 'H 20 25400 001' and 'HE' status. It includes a 'Hematoxilin Eosin' dropdown and 'Request' and 'Delete' buttons.

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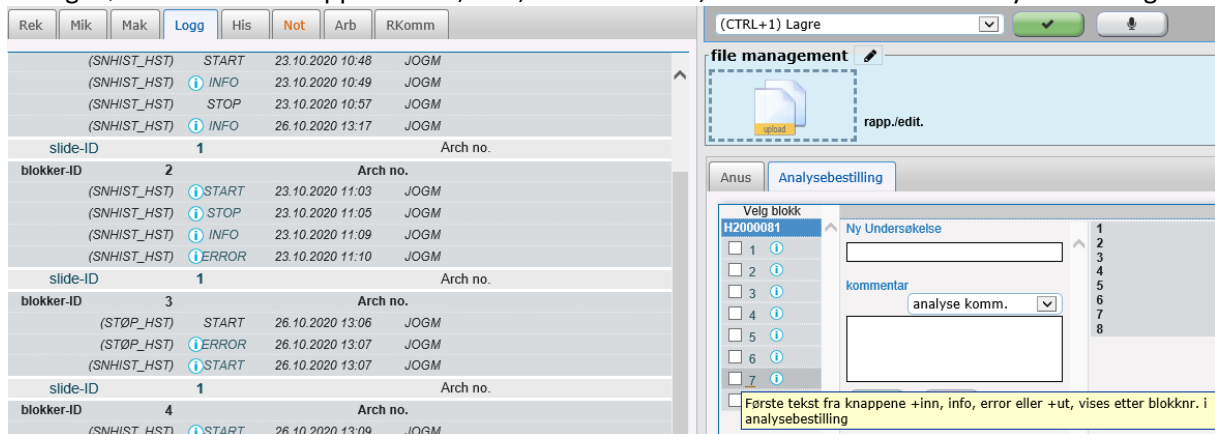
The screenshot shows the UniApp interface with a log of events on the left and a file management window on the right. The log shows several events including START, INFO, STOP, and ERROR for different blocks. The file management window shows a list of blocks and a selected block with a comment field.

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The screenshot shows the UniApp interface with a control panel on the left and a history log on the right. The control panel includes buttons for 'inn', 'info', 'Error', and 'ut', along with a status bar showing 'STØP (STØP\_HST) missing START/STOP'. The history log shows a list of events with columns for Date, Time, Process Point, Workplace, User, and Device.

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